

USING TIME CONTROLLER FOR FLUORESCENCE MICROSCOPE

Hongwei Cheng 8/22/01, commented by TCH 1/27/02

How to use it:

1. Press "ON/OFF" key to turn the timer on.
2. Press "CD" key to start CD program. The timer will automatically turned off itself 15 minutes later. In turn, the UV supply unit will be powered off.

Notes:

1. You are not supposed to use Time Controller every time. You need it just in the case that your looking time is less than 15 minutes and you don't want to wait until 15 minutes to turn it off.
2. If your looking time is more than 15 minutes, what you need to do is just to press "ON/OFF" key to turn the timer off.
3. Due to the 15 minutes turn on and turn off interval rule, we always set up the CD program on this point. So usually, you don't need to set it up again.

If you have any problems in operating the timer, please see Hongwei or TCH